

SRR&CVR GOVT. DEGREE COLLEGE(A), VIJAYAWADA

Department of Biochemistry

Hands on Training

Activity: Guest lecture and hands on training by Mrs. Dorka Vijaya, Lecturer in Biochemistry, GDC Women, Guntur.

Date: 13/03/2019

Description: The department Biochemistry, SRR&CVR GDC, Conducted hands on training on protein separation by SDS PAGE. Mrs. Dorka Vijaya from GDC Women, Guntur acted as resource person, demonstrated the technique, and separated the protein "serum albumin". Students also practiced the gel loading and sample loading techniques by hands on training. We thank our principal Dr. Joshi sir for giving permission to buy the SDS PAGE kit and to conduct this hands on training programme. Faculty from department of zoology and microbiology also participated in this training programme.



SRR&CVR GOVT. DEGREE COLLEGE(A), VIJAYAWADA
Department of Biochemistry



SRR&CVR GOVT. DEGREE COLLEGE(A), VIJAYAWADA
Department of Biochemistry



SRR&CVR GOVT. DEGREE COLLEGE(A), VIJAYAWADA

Department of Biochemistry



SRR & CVR GOVT. DEGREE COLLEGE

(Autonomous)

NAAC accredited with 'B' Grade

Machavaram, Vijayawada – 520 004, Krishna District.

Cell: 94922 34488 Ph: 0866-2430060 Fax: 0866-2441092 www.srrcvt.org srrandcvt@gmail.com



Dr. Velaga Joshi, Principal

M.A.(PhD), M.A. (Eng), M.A. (MLC), LL.B., M.Phil., Ph.D.

Dt: 12-03-2019.

To,
Dorka Vijaya Kumari.B
Assistant Professor,
Department of Biochemistry,
GCW (Autonomous),
Guntur.

Respected Sir,

Sub: Invitation to act as a resource person to demonstrate & explain the technique SDS PAGE in Department of Biochemistry at S.R.R& C.V.R Govt (A) DEGREE COLLEGE.

The Department of Biochemistry is proposing to conduct workshop on 13th March 2019 to motivate and to create awareness among under graduate students on the technique of Gel electrophoresis (SDS PAGE).

Your expertise and experience in this technique will be an excellent resource to our students and it is honor and pleasure to invite you to demonstrate the technique.

On behalf of our life sciences faculty I am requesting you to kindly accept our invitation and I look forward for your positive response.

Thanking You,

Yours sincerely,

PRINCIPAL
SRR & CVR GOVT. DEGREE COLLEGE
(Autonomous)
Machavaram, VIJAYAWADA-520 004.

Laemmli SDS PAGE

Fanglian He

Carnegie Institution at Stanford

[Abstract] Sodium dodecyl sulfate polyacrylamide gel electrophoresis (SDS-PAGE) is used to separate proteins with relative molecular mass no smaller than 10 KD. Very small proteins (<10 KD) are difficult to resolve due to low ability of binding to SDS, which can be solved by gradient gels or using different electrophoresis conditions, like Tricine-SDS-page. The basic Laemmli SDS PAGE procedure is described here.

Materials and Reagents

1. Pre-stain Protein MW marker (Bio-Rad Laboratories)
2. TEMED (Life Technologies, Gibco[®])
3. Ammonium persulfate (Sigma-Aldrich)
4. SDS (Research Organics INC)
5. 30% Acrylamide stock (37.5: 1 acrylamide: bisacrylamide) (Bio-Rad Laboratories)
6. Bromophenol Blue (Thermo Fisher Scientific)
7. Tris Base (Calbiochem-Behring)
8. Glycine (EM Science)
9. EDTA (Amresco)
10. Glycerol (EM Science)
11. Isopropanol
12. Tris-HCl (pH 6.8)
13. β -mercaptoethanol (Sigma-Aldrich)
14. 10x running buffer (see Recipes)
15. 2x SDS protein sample buffer (see Recipes)

Equipment

1. Protein mini gel cassettes (Bio-Rad)
2. Heating block module
3. Table-top centrifuge
4. Power supply
5. Gloves

6. Filter paper

Procedure

A. Making SDS-PAGE gel

1. Clean and completely dry glass plates, combs, and spacers are required.
2. Assemble gel cassette by following manufacturer instructions.
3. Prepare 10% lower gel (separating gel) by adding the following solutions (wear gloves when prepare gel solution) (total volume= 5 ml)
2 ml ddH₂O
1.67 ml 30% acrylamide/Bis
1.25 ml 1.5 M Tris (pH 8.8)
25 µl 20% SDS
25 µl 10% ammonium persulfate (make it fresh and store at 4 °C up to a month)
2.5 µl TEMED (add it right before pour the gel)
Note: Change ratio of ddH₂O to 30% acrylamide/Bis to get different percentage of separating gel.
4. To avoid polymerization, after adding TEMED, mix well and quickly transfer the gel solution by using 1 ml pipette to the casting chamber between the glass plates and fill up to about 0.7 cm below the bottom of comb when the comb is in place.
5. Add a small layer of isopropanol to the top of the gel prior to polymerization to straighten the level of the gel.
6. Once the gel has polymerized, start to prepare stacking gel (5%) by adding the following solutions (total volume= 3 ml)
2.068 ml dH₂O
0.506 ml 30% acrylamide/Bis
0.375 ml 1 M Tris (pH 6.8)
15 µl 20% (w/v) SDS
15 µl 10% ammonium persulfate
1.5 µl TEMED (add it right before the gel is poured)
7. Remove the isopropanol layer by using filter paper. Rinse the top layer of the gel with ddH₂O and dry off as much of the water as possible by using filter paper.
8. Add TEMED and mix the stacking gel solution well. Quickly transfer the gel solution by using a 1 ml pipette till the space is full, and then insert the appropriate comb.
9. Allow the top portion to solidify and then carefully remove the comb.

Note: The gels can be stored with the combs in place tightly wrapped in plastic wrap and put in a second container with wet tissue towel (keep the gels moist) at 4 °C for 1 to 2 weeks.

B. Sample preparation

1. Prepare same amount of protein samples according to BCA assay result, see [BCA \(bicinchoninic acid\) protein assay](#).
2. Add the same volume of 2x protein sample buffer to each protein sample, mix and boil the samples at 95 °C heating block module for 10 min.
3. Spin the samples at the maximal speed for 1 min (samples from some tissue/cell sources may need longer spin) in tabletop centrifuge and leave the samples at room temperature until you are ready to load onto the gel.

Note: Can store extracted protein samples (containing sample buffer) at -20 °C and reheat at 95 °C for 5 min when used the following time.

C. Electrophoresis

1. Remove the gel cassette from the casting stand and place it in the electrode assembly with the short plate on the inside. Press down on the electrode assembly while clamping the frame to secure the electrode assembly and put the clamping frame into the electrophoresis tank.
2. Pour some 1x electrophoresis running buffer into the opening of the casting frame between the gel cassettes. Add enough buffer to fill the wells of the gel. Fill the region outside of the frame with 1x running buffer.
3. Slowly load the same amount of protein samples into each well as well as load 10 µl of protein MW marker.
4. Connect the electrophoresis tank to the power supply.

D. Protein detection

If protein of interest is about 0.2 µg or more in the sample, typically use Coomassie blue staining (see [Coomassie blue staining](#)). Otherwise, use silver staining (silver staining), which is more sensitive and can detect as little as 5 ng protein.

Recipes

1. 10x running buffer
30.3 g Tris-base
144.0 g glycine

10.0 g SDS

Completely dissolve in about 800 ml ddH₂O and then more ddH₂O up to 1 liter.

2. 2x SDS protein sample buffer

1.25 ml 1 M Tris-HCl (pH 6.8)

4.0 ml 10% (w/v) SDS

2.0 ml glycerol

0.5 ml 0.5 M EDTA

4 mg bromophenol blue

0.2 ml β-mercaptoethanol (14.3 M)

Bring the volume to 10 ml with ddH₂O.

References

1. Laemmli, U. K. (1970). [Cleavage of structural proteins during the assembly of the head of bacteriophage T4.](#) *Nature* 227(5259): 680-685.

PROTEIN SEPERATION

BY SDS PAGE

13/3/19

70

The dept. of Biochemistry has conducted workshop on SDS PAGE Technique on 13/3/19 for this Smt. DRCas Vijaya, Assistant professor, Dept. of Biochemistry, Guntur women's degree college acted as a resource person, and performed the technique, seperated Serum sample & Ladder proteins sample from the dept. of Zoology. Smt. G. Vani garu, Smt. Parveen Madan garu from botany dept. Smt. D. Jyothi garu participated and all 1st, 2nd, 3rd Year students of MBC group were ~~also~~ engaged in the workshop. 13/3/19.

Smt. D. Jyothi
I/c biochemistry

Smt. D. Jyothi

G. Parvathi

Smt. G. Vani

Smt. Parveen Madan

G. Nagalakshmi

IIIrd B.Sc (MBC)

Nagalakshmi

M. Ratnakumari

IIIrd B.Sc [MBC]

M. Ratnakumari

S. Kunti

IIIrd B.Sc (MBC)

S. Kunti

B.S. Ketharin

IIIrd B.Sc (MBC)

B.S. Ketharin

A. Poovanchander

IIIrd B.Sc (MBC)

A. Poovanchander

Student Signature

71

1. M. Chaitanya Hanish	III rd BSc (MBC)	(Signature)
2. R. Durga Rajeswari	" "	R. D. Rajeswari
3. B. Teja Naga Sai	I st BSc (MBC)	B. Teja Naga Sai
4. G. Surya Prakash	I st B.Sc (MBC)	G. Surya Prakash
5. G. Neeraja	I st B.Sc (MBC)	G. Neeraja
6. K. Parvathi	I st B.Sc (MBC)	K. Parvathi
7. L. Lalitha Bhargavi	I st B.Sc (MBC)	L. Lalitha Bhargavi
8. A. Nazma	III BSc (MBC)	A. Nazma
9. M. Vyshnavi	III BSc (MBC)	M. Vyshnavi
10. A. Raghunadh Reddy	I BSc (MBC)	A. Raghunadh Reddy
11. T. Yasaswini Sri Sai	III BSc (MBC)	T. Yasaswini Sri Sai
12. K. Vijaya Krishna	III BSc (MBC)	K. Vijaya Krishna
13. K. Bhavani	III BSc (MBC)	K. Bhavani
14. MD. Shahid Ali	III BSc (MBC)	MD. Shahid Ali